# FEATHER DISPOSABLE MICROTOME BLADES & HOLDER No.240D

- FEATHER MIKROTOMKLINGEN & HALTER No.240D
- LAMES POUR MICROTOME ET PORTE-LAME FEATHER No. 240D
- CUCHILLAS PARA MICRÓTOMO Y SOPORTE FEATHER No. 240D
- LAME PER MICROTOMO E PORTALAMA FEATHER No. 240D
- LÂMINAS PARA MICRÓTOMO E SUPORTE FEATHER N.º 240D
- FEATHER MICROTOOMMESJES & HOUDER Nr. 240D

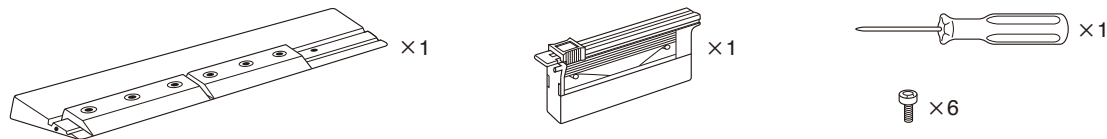
## INSTRUCTIONS FOR USE

- GEBRAUCHSANWEISUNG
- INSTRUCTIONS D'UTILISATION
- INSTRUCCIONES DE USO
- ISTRUZIONI PER L'USO
- INSTRUÇÕES DE UTILIZAÇÃO
- GEBRUIKSAANWIJZING



## CONTENTS

- INHALT
- CONTENU
- CONTENIDOS
- CONTENUTI
- CONTEÚDO
- INHOUD

## HOW TO CLAMP THE DISPOSABLE MICROTOME BLADE

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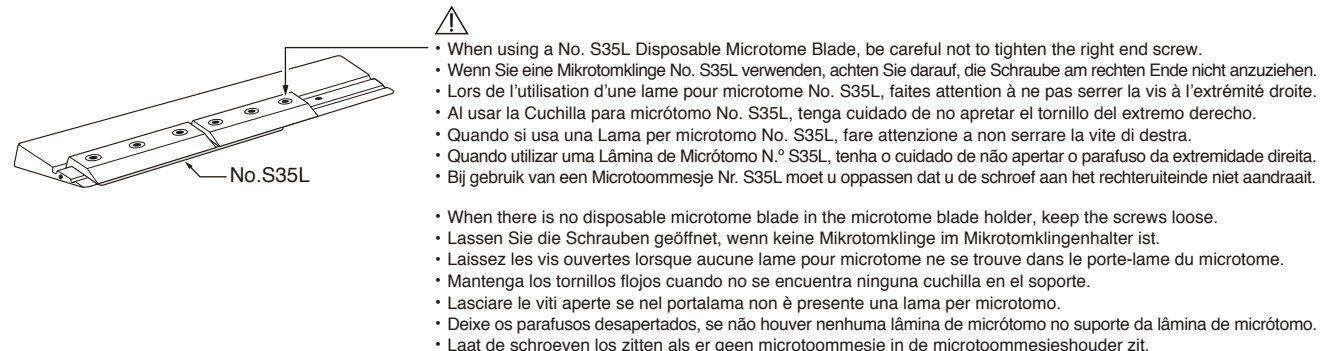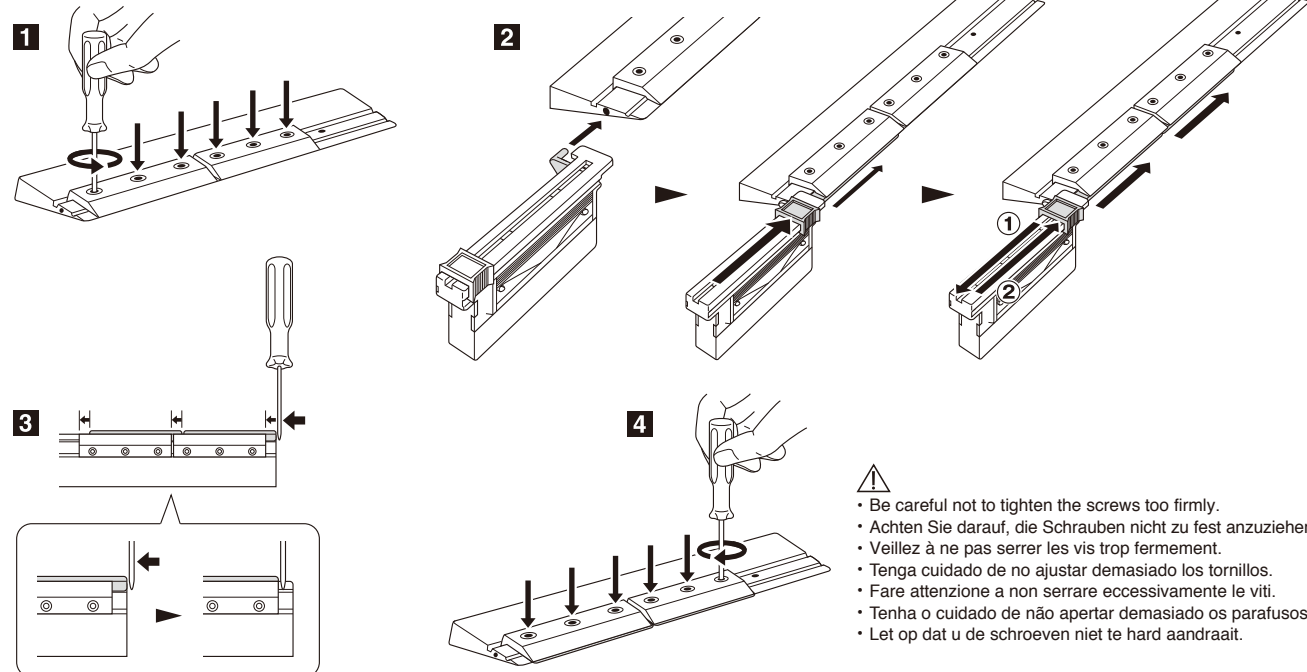
## CAUTION
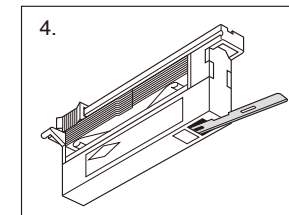
### PRECAUTIONS FOR USE

1. The product is only for medical use. Disposable microtome blades are extremely sharp, so please use them with care. Used microtome blades can be safely kept in the part of the dispenser marked accordingly.
2. To insert a new disposable microtome blade, please use the insertion device on the microtome blade holder or dispenser.
3. Make sure to fix the microtome blade holder at a clearance angle of at least 5°.
4. Used microtome blades can be safely kept in the part of the dispenser marked accordingly.
5. The disposable microtome blades have been treated with anti-corrosion oil. This will not cause any problems for the sectioning process, so removal of the oil is not necessary.
6. Disposable microtome blades will quickly become dull if used for paraffin blocks which have not been well infiltrated with paraffin, for sectioning non-decalcified tissues or tissues having minute structures.

### SAFETY AND CARE INSTRUCTIONS

1. Wipe the surface of the microtome blade holder daily using machine oil. Although the holder is made of alloyed tool steel and is rust-resistant, moisture and fingermarks could cause rust.
2. Handle with care in order not to damage the precision ground edge of the microtome blade holder.
3. Prior to storing the microtome blade holder for a longer period of time, fully disassemble all the parts and apply some oil to all surfaces, including the grooves, fixing plates and screw holes. **DE**

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